

Instructions Isohelix Stabilising and Lysis kits -- DSK-3/DSK-50

STORE COMPONENTS AT CORRECT TEMPERATURES ON ARRIVAL

Product Details.

April 2007

The Isohelix Stabilising and Lysis kits are designed to fully stabilise your DNA (taken from swabs and other source material) from enzymatic activity that occurs naturally after Buccal sampling. The kit will store the samples at room temperature and shows no loss of stability even after 7 months of data collected to date. The DNA LS Kit also contains a component to fully lyse the cells as well as enzymatic inhibitors, releasing all the DNA and maintaining the structural integrity.

- * Stabilises your DNA
- * Room Temperature storage
- * Long term storage for DNA

- * Improves DNA yield and Purity
- * Lyses the cells
- * Pre- treatment prior isolation

Storage and Stability

The DNA Isolation kits are shipped at ambient temperature, on arrival **PLEASE NOTE** that some of the components should be stored according to the table below upon arrival. The kit is stable for 2 Years if stored as instructed.

Vial	Storage Condition
Solution LS	Room Temperature Approx. +20°C to +22°C
Solution PK	+4°C

Kit Contents.

Catalogue No: DSK-3 - Trial kit

Vial	No.	Qty	Contents
Solution LS	1 x	1.5 ml	Lysis/Storage Buffer
Solution PK	1 x	100 µl	Proteinase K buffer

One kit contains enough material for 3 extractions.

Catalogue No: DSK-50

Vial	No.	Qty	Contents
Solution LS	1 x	25 ml	Lysis/Storage Buffer
Solution PK	1 x	1 ml	Proteinase K buffer

One kit contains enough material for 50 extractions.

Before you start:

If a precipitate has formed in Solution LS incubate the tube/bottle at 60°C for few minutes.

DNA stabilising and Lysis from the swab

1. Add 500 µl of LS solution to the tube containing the swab.
2. Add 20 µl of PK. Solution, Vortex briefly

NB: The DNA can now be extracted or the sample can be kept for at least 7 months at room temperature.

Other Products for DNA sample handling

Isohelix Buccal Swabs:

High yields, blood alternative, reproducible, easy to use, different formats for various extraction methods

Isohelix DNA Isolation Kits

DDK 3/50

Optimised for buccal cells, no organic solvents, reduced steps, HT formats, No columns or filters.

Isohelix DNA quality check

DSK/50

PCR kit to confirm quality of DNA prior onward experimentation.

Safety and Use of the DSK kits

The DSK kits are intended for use by qualified professionals trained in potential laboratory hazards and good laboratory practise. If direct information is not available on any of our compounds this should not be interpreted as an indication of product safety.

This kit has been designed solely for research use only

Isohelix is a division of Cell Projects Ltd

For Swab and DNA isolation queries info@isohelix.com - www.isohelix.com
For Molecular Biology Solutions www.cellprojects.com